



FitMe™ Lab Jackets and Coats Care Instructions

Autoclave:

1. Using only one garment at a time, fold loosely and place into the autoclave
2. Set autoclave on 180 degrees
3. Remove garment from autoclave before drying cycle
4. Shake to remove wrinkles
5. Hang to dry
6. WARNING: DO NOT CHEMICLAVE

Laundry:

1. Machine wash in warm water with bleach
2. Shake to remove wrinkles
3. Hang to dry
4. WARNING: DO NOT USE A COMMERCIAL DRYER

Please Note:

- Dukal does not guarantee that either autoclave or laundry will kill specific viruses or bacteria
- Autoclave machines must be maintained and calibrated regularly
- FitMe™ Lab Coats and Jackets can be autoclaved or washed up to 5 times
- If you notice any tears, particularly at the seam, discard the garment



Find our full line of products at [dukal.com](https://www.dukal.com)

Dukal, LLC
2 Fleetwood Ct, Ronkonkoma, NY 11779

Phone: 1 631 656 3800
Toll Free: 1 800 243 0741

Email: customerservice@dukal.com
To Place Orders: orders@dukal.com